### Alaska Scientific Crime Detection Laboratory

#### DataMaster Check In Form

Issued: 1/13/2017 Version: DMCIF 2017 R0 Effective: 1/17/2017 Status: Active Instrument S/N /00667 Analyst Derch Watton 02/25/20 Date 1. Reason for instrument return: 3W 02/25/20 Returned Fo to factory for refurbish and because it would not recall lamp and cooler voltages when shut down. 2. Change instrument and case location in JusticeTrax Breath Alcohol Database 3. Ensure change in instrument status form is in JusticeTrax Not applicable DW 02/26/20 4. Upload 5. Update agency information and polling status in DM Host Calling Program 6. Check for VOCs

#### From Intoximeters

7. Intoximeters Repair Paperwork (attached)

Instrument Sent to Intoximeters on 01/13/20 for repair.
Repair completed on 02/13/20 and shipped back to lab on 02/17/20. Leceived by lab on 02/19/20. JM 03/03/20

# Intoximeters

Experience · Service · Integrity

### Service Record

SRO Number: SRO-038483

**Received Date:** 01/15/2020

Completion Date: 02/13/2020

SRO Type: REPAIR

SRO Description: DMT Repair

**Customer ID:** C000003651

Customer PO: 20200108DJW

Contact: Derek Walton

Email: derek.walton@alaska.gov

Ship To:

Units on SRO

Alaska Scientific Crime Detection Laboratory

4805 Dr MLK Jr Ave Anchorage AK 99507

USA

LEGDMT LEGACY DMT

Service Line:	**All Instruments are Calibrated to Factory Specifications**					
Unit Type:	(LEGDMT) LEGACY DMT					
Serial Number:	50100667					
Reason:	DMT Functional					
	No Boot					
Reason Notes:	: Customer Reported Issue: Will not recall correct lamp and cooler voltages after restart, controller board with jumper may be needed.					
	TEC: See SRO notes prior to repairing.					
Resolution:	Prev M					
	planned					
	Installed TB-30 modification Leak checked dry gas regulator Rebuilt filter assembly and added epoxy to motor wires Optimized detector Cleaned and adjusted 5-way valve Adjusted tech screen settings and RFI sensitivity Added o-ring to end of breath tube Replaced leaking 'Vapor In' panel mount connector					

Parts:	Qty	Part Number	Part Description
	1	001199-00	DMT ELEC REGULATOR V LDO 3.3V OUT 30V IN
	1	59-0042-00	DMT HRDW CPC PANEL MNT MALE 26002
	1	90-8000-00	CONTRACT PREV MAINT DMT

# Order Packing Slip Report

2/17/2020 2:47:19 PM

Intoximeters

Experience · Service · Integrity

Packing Slip: 198319

Pack Date: 2/17/2020

SRO: SRO-038483

From: Intoximeters, Inc

2081 Craig Road St Louis MO 63146

USA

Customer: C000003651

Ship To: 1

Alaska Scientific Crime Detection

Alaska Department of Public Safety

4805 Dr MLK Jr Ave Anchorage AK 99507

USA

Laboratory 4805 Dr MLK Jr Ave

Anchorage AK 99507

USA

Customer PO: 20200108DJW

Ship Via: FDX

FED EX

Weight: 37.00

Delivery Terms: 2D

FOB: ORIGIN

# Packages: 1

**Customer Ship Account:** 

Line Number	Item		dered antify	Shipment Quantity	Backorder Quantity				
1	LEGDMT	LEGACY DM	Т		50100667				
1	001199-00	EA		1.00	0.00				
DMT ELEC REGULATOR V LDO 3.3V OUT 30V IN									
2	59-0042-00	EA		1.00	0.00				
DMT HRDW CPC PANEL MNT MALE 26002									
3	90-8000-00	EA		1.00	0.00				
CONTRACT PREV MAINT DMT									