Alaska Scientific Crime Detection Laboratory

DataMaster Check In Form

	ssued: 1/13/2017 ffective: 1/17/2017	Version: DMCIF 2017 R0 Status: Active
Instr	ument S/N <u>/104/3</u>	Received From Internation
Date_	04/22/21	Analyst Deren Walton
1.	Reason for instrument return: Returnished by formulator the	etory repaired for tanh interference and neont status messages
2.	Change instrument and case location in	n JusticeTrax Breath Alcohol Database
3.	Ensure change in instrument status form	
4.	Upload	Instrument was not in the field
5.	Update agency information and polling	
6.	Check for VOCs	
	From Intoximeters	
7.	Intoximeters Repair Paperwork (attache	d) 🗹

Experience - Service - Integrity

Service Record

SRO Number: SRO-045230

Received Date: 03/30/2021

Completion Date: 04/09/2021

SRO Type: REPAIR

SRO Description: DMT Preventive Maintenance

Customer ID: C000003651

Customer PO: 033021-Walton

Contact: Derek Walton

Email: derek.walton@alaska.gov

Ship To:

Units on SRO

1	LEGDMT	LEGACY DMT	
	1	1 LEGDMT	1 LEGDMT LEGACY DMT

Service Line:	**All Instruments are Calibrated to Factory Specifications**
Unit Type:	(LEGDMT) LEGACY DMT
Serial Number:	50100413
Reason:	DMT Functional
	Lock Up
Reason Notes:	Customer Reported Issue: Tank interference, simulator timeout using dry gas tank
Resolution:	Adjusted
	Elec adj
	Adjusted Tech screen settings Replaced filter wheel solenoid mounting screws Replaced stepper motor adjustment lock down screw Replaced filter wheel motor pinion Cleaned, lubricated, & adjusted filter wheel moving parts Optimized detector position

Parts:	Qty	Part Number	Part Description				
	1	001199-00	DMT ELEC REGU	JLATOR V LDO 3.3V OUT 30V IN			
Reason:	n: Preventive Maintenance						
	planned						
Reason Notes:	Preventative Maintenance contract						
Resolution:	Prev M						
	planned						
Resolution Notes:	Tightened dry gas port panel mount nut Replaced dried/cracking breath tube inlet o-ring Trimmed .5 inches of unheated tubing Adjusted tech screen settings Adjusted RFI sensitivity Sealed mirror blocks Installed Torx screws on sample chamber Installed TB-30 DMT Digital Pots Mod Adjusted 5 way valve solenoid position Tested modem connectivity Adjusted air volume Tested and verified coin battery: 3.147vdc Parts: 1-001199-00 Voltage Reg						
Cal. solution (0.100) Lot#21080 Bot#1206 exp.03/08/2023; Dry gas: AG919705, Tank 029, Exp. 16JUL2022; AG004303, Tank 053, Exp. 12Feb2022							
	1	90-8000-00	CONTRACT PRE	CONTRACT PREV MAINT DMT			

Order Packing Slip Report

Intoximeters
Experience · Service · Integrity

Pack Date: 4/14/2021

4/14/2021 1:29:21 PM

Packing Slip: 225918

SRO: SRO-045230

From: Intoximeters, Inc

2081 Craig Road St Louis MO 63146

USA

Customer: C000003651

Alaska Department of Public Safety

4805 Dr MLK Jr Ave Anchorage AK 99507

USA

Ship To: 1

Alaska Scientific Crime Detection

Laboratory

4805 Dr MLK Jr Ave Anchorage AK 99507

USA

Customer PO: 033021-Walton

Ship Via: FDX

FED EX

Weight: 37.00

Delivery Terms: 2D

FOB: ORIGIN

Packages: 1

Customer Ship Account:

Line Numl	oer Item	UM	Ordered Quantity	Shipment Quantity	Backorder Quantity		
1	LEGDMT	LEGACY	Y DMT		50100413		
1	001199-00	EA		1.00	0.00		
DMT ELEC	C REGULATOR	V LDO 3.3	V OUT 30V IN				
2	90-8000-00	EA		1.00	0.00	iluò	
2 90-8000-00 EA 1.00 0.00 juid CONTRACT PREV MAINT DMT ONIGHT							
						DM	