Alaska Scientific Crime Detection Laboratory

DataMaster Check In Form

	D 0.001, 10.000.	O
Issued: 1/13/2017		Version
Effective: 1/17/2017		

ersion: DMCIF 2017 R0/ Status: Active

Instrument S/N	Received From Into Kineters
Date02/14/20	Analyst Derew Walton

1. Reason for instrument return:

Instrument repaired	by Intoximeters for	bad 5 way valve and
unably to pass line	easty	V
	/	

- 2. Change instrument and case location in JusticeTrax Breath Alcohol Database
- 3. Ensure change in instrument status form is in Justice Trax [Not appliable) WOZ/14/20
- 4. Upload | Not applicable DW 02/14/20
- 5. Update agency information and polling status in DM Host Calling Program \[\int_{Not} appliable \) W 02/14/20
- 6. Check for VOCs 1 Not applicable 02/14/20

From Intoximeters

7. Intoximeters Repair Paperwork (attached)

Instrument sent to Intoxincters on 01/13/2020 and received by Intoximeters on 01/15/2020. Instrument returned from Intoximeters on 02/14/2020 (date of checkin). DM 03/02/2020

Intoximeters

Experience · Service · Integrity

Service Record

SRO Number: SRO-038485

Received Date: 01/15/2020

Completion Date: 01/28/2020

SRO Type: REPAIR

SRO Description: DMT Repair

Customer ID: C 000003651

Customer PO: WTY03

Contact: Derek Walton

Email: derek.walton@alaska.gov

Ship To:

Units on SRO

Alaska Scientific Crime Detection Laboratory 4805 Dr MLK Jr Ave Anchorage AK 99507

LEGDMT

LEGACY DMT

USA

Service Line:	**All Instruments are Calibrated to Factory Specifications**		
Unit Type:	(LEGDMT) LEGACY DMT		
Serial Number:	50100367		
Reason:	DMT Functional		
	Erratic		
	Customer Reported Issue: did not pass linearity when received, blank errors and adjustment errors on attempted calibration, 5 way valve also sticking, replaced 5 way valve block assembly with new parts received from Intoximeters - gets pump errors with simulator use preventing calibration - flow voltage when sim valve is closed is too low, suspect issue with new 5 way valve block assemblies purchased from Intoximeters.		
Resolution:			
4.1.21.424 1.1.1.1.1.1.1.1.1.1.1.1.1.1.1.1.1.1.1.			
	Adjusted touch screen Replaced 5-way valve Cleaned filter optics Adjustment and accuracy tested Adjusted RFI sensitivity		

Parts:	Oty	Part Number	Part Description
	1	59-0018-00	DMT HRDW ASSY VALVE 5WAY 12VDC 41505

Order Packing Slip Report

1/31/2020 7:47:09 AM

Intoximeters

Pack Date: 1/31/2020

Experience · Service · Integrity

1,01,2020 /.....09 /...

Packing Slip: 196829

SRO: SRO-038485

From: Intoximeters, Inc

2081 Craig Road St Louis MO 63146

USA

Customer: C000003651 Ship To: 1

Alaska Department of Public Safety

4805 Dr MLK Jr Ave Anchorage AK 99507

USA

Alaska Scientific Crime Detection

Laboratory

4805 Dr MLK Jr Ave Anchorage AK 99507

USA

Customer PO: WTY03

Ship Via: FDX

FED EX

Weight: 38.00

Delivery Terms: 2D

FOB: DESTINATION

Packages: 1

Customer Ship Account:

Line Number	fem	UM Ordered Quantity	Shipment Quantity	Backorder Quantity	
1	LEGDMT	LEGACY DMT		50100367	
1	59-0018-00	EA	1.00	0.00	lec'd 02/14/20 DM
DMT HRDW A	SSY VALVE	5WAY 12VDC 41505			peca 0=1. 100 m