

Alaska Scientific Crime Detection Laboratory

DataMaster Check In Form

Issued: 1/13/2017
Effective: 1/17/2017

Version: DMCIF 2017 R0
Status: Active

Instrument S/N 100 678

Received From Intoximeter

Date 09/15/21

Analyst Derek Watten

1. Reason for instrument return:

Rebuilt by manufacturer, repaired for crashing on startup

2. Change instrument and case location in JusticeTrax Breath Alcohol Database

3. Ensure change in instrument status form is in JusticeTrax

4. Upload

*NA Instrument was not in the field
DMW 09/15/21*

5. Update agency information and polling status in DM Host Calling Program

6. Check for VOCs

From Intoximeters

7. Intoximeters Repair Paperwork (attached)

SRO Number: SRO-047143

Received Date: 07/26/2021

Completion Date: 08/18/2021

SRO Type: REPAIR

SRO Description: DMT Repair/Preventative Maintenance

Customer ID: C000003651

Customer PO: 07262021-Lee

Contact: Tatiana Lee

Email: tatiana.lee@alaska.gov

Ship To:

Alaska Department of Public Safety
Procurement & Supply Section
4805 Dr. MLK Jr. Avenue
Anchorage AK 99507

Units on SRO

1	LEGDMT	LEGACY DMT
---	--------	------------

Service Line: 1	**All Instruments are Calibrated to Factory Specifications**
-----------------	---

Unit Type:	(LEGDMT) LEGACY DMT
Serial Number:	50100678
Reason:	DMT Functional
	No Boot
Reason Notes:	Customer Reported Issue: Crashes on Startup
Resolution:	Update
	Update
Resolution Notes:	Found TicketDB file corrupted. Downloaded files off of SD Card and reformatted Memory Card. Loaded files back onto SD Card minus the corrupted TicketDB file. Instrument boots up correctly.

Parts:	Qty	Part Number	Part Description
	1	90-8000-00	CONTRACT PREV MAINT DMT
Reason:	Preventive Maintenance		
	planned		
Reason Notes:	Maintenance of Filter Wheel Assembly, Sample Chamber, and 5-Way Valve.		
Resolution:	Prev M		
	planned		
Resolution Notes:	Removed filter wheel assembly for cleaning and lubrication. Cleaned & adjusted filter wheel solenoid. Updated filter wheel solenoid mounting screws. Updated stepper motor adjustment lock down screw. Replaced filter wheel motor pinion. Cleaned, lubricated, & adjusted filter wheel moving parts. Removed, inspected, and applied metallic seal tape to mirror blocks of Sample Chamber. Replaced discolored tubing and inspected 5-Way Valve. Verified Modem with laptop and DMT Tech Software using phone cord.		
	1	001199-00	DMT ELEC REGULATOR V LDO 3.3V OUT 30V IN
Reason:	DMT Functional		
	DigPots		
Reason Notes:	Requires Digital Pots Mod.		
Resolution:	Update		
	DMTVR3		
Resolution Notes:	Removed Control Board and installed TB-30 DMT Digital Pots Mod. Optimized detector position. Adjusted Tech Screen settings.		

Order Packing Slip Report

Intoximeters
Experience · Service · Integrity

8/26/2021 9:23:46 AM

Packing Slip: 235294

Pack Date: 8/26/2021

SRO: SRO-047143



From: Intoximeters, Inc
2081 Craig Road
St Louis MO 63146
USA

Customer: C000003651

Ship To: 2

Alaska Department of Public Safety
4805 Dr MLK Jr Ave
Anchorage AK 99507
USA

Alaska Department of Public Safety
Procurement & Supply Section
4805 Dr. MLK Jr. Avenue
Anchorage AK 99507
USA

Customer PO: 07262021-Lee

Ship Via: FDX

FED EX

Weight: 38.00

Delivery Terms : 2D

Shipping Terms : ORIGIN

Packages: 1

Customer Ship Account :

Line Number	Item	UM	Ordered Quantity	Shipment Quantity	Backorder Quantity
1	LEGDMT	LEGACY DMT			50100678
1	90-8000-00	EA		1.00	0.00
	CONTRACT PREV MAINT DMT				
2	001199-00	EA		1.00	0.00
	DMT ELEC REGULATOR V LDO 3.3V OUT 30V IN				